



AQUA-TECH MEDIA PREP

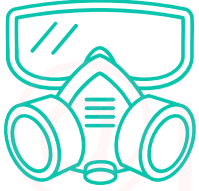

WEAR APPROPRIATE DUST MASK, GOGGLES, GLOVES, AND PROTECTIVE CLOTHING AT ALL TIMES.

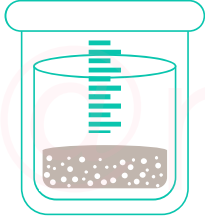

**PREPARE MEDIA IN A THIN LAYER
SPREAD EVENLY ACROSS THE BOTTOM OF A VERY
DEEP STAINLESS STEEL OR GLASS CONTAINER**



**USE A VACUUM OVEN, JACKETED REACTOR OR
OTHER AMBIENT HEAT SOURCE.**

**DO NOT USE CONVECTION OVENS OR DEHYDRATORS
WITH FANS.**

***SERIOUSLY, DO NOT USE MOVING AIR TO BAKE YOUR MEDIA!**



BAKE 200C FOR 6-8 HOURS



**WEIGH YOUR DRY MEDIA INDIVIDUALLY, THEN
COMBINE THEM IN A LARGE, AIRTIGHT CONTAINER
NO MORE THAN 1/3 FULL.**



**SIFT AND STIR THE DRY MEDIA TO BREAK UP ANY
CHUNKS WHILE TOTALLY BLENDING THE MIXTURE.**



**RECORD THE PREPARATION DATE, CHEMICAL
CONTENT - INCLUDING BATCH NUMBERS AND
WEIGHTS - ON THE EXTERIOR OF THE CONTAINER, AS
WELL AS IN A PERMANENT LOG FOR FUTURE
REFERENCE.**